


SONY

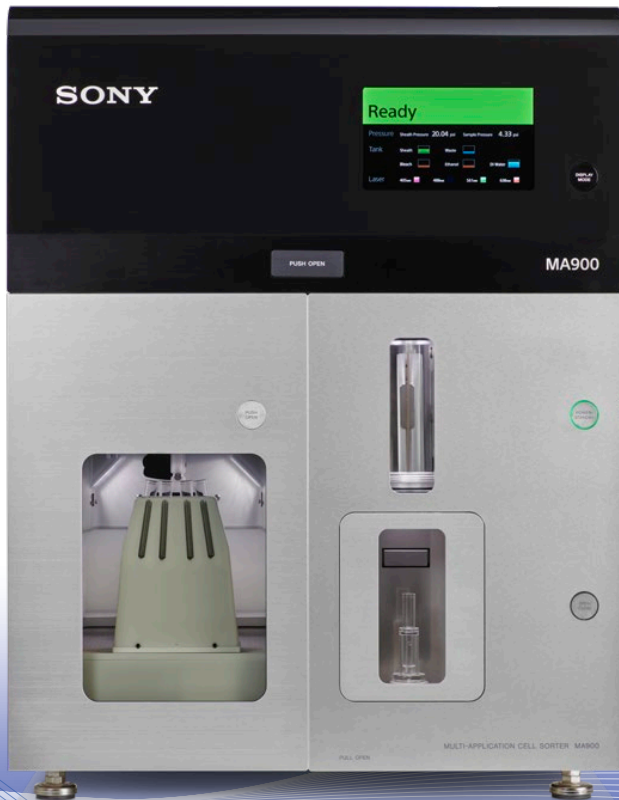

MA900 Multi-Application Cell Sorter

Sony Biotechnology Inc.

Fluorochrome and Filter Guide

Fluorochrome		FL1 525/50	FL2 585/30	FL3 617/30	FL4 695/50	FL5 785/60	
EGFP		●					
FITC		●					
Alexa Fluor® 488		●					
EYFP		●					
mCitrine		●					
CFSE		●					
PE			●				
dsRed			●				
tdTomato			●				
mCherry				●			
PE-Texas Red®				●			
PE/Dazzle 594				●			
Propidium iodide				●			
mPlum					●		
7-AAD					●		
PE-Cy™5					●		
PerCP					●		
PE-Cy5.5					●		
PerCP-Cy5.5					●		
PerCP-eFluor 710					●		
PE-Cy7						●	
	FL6 450/50	FL7 525/50	FL8 585/30	FL9 617/30	FL10 665/30	FL11 720/60	FL12 785/60
BD Horizon Brilliant Violet (BV421)	●						
Alexa Fluor® 405	●						
DAPI	●						
Pacific Blue™	●						
mCFP	●						
Hoechst 33342	●						
AmCyan		●					
BV510		●					
BV570			●				
BV605				●			
BV650					●		
BV711						●	
BV785							●
APC					●		
Alexa Fluor® 647					●		
APC-Cy5.5						●	
Alexa Fluor® 700						●	
APC-Cy7							●
APC-Alexa Fluor® 750							●

Filter Guide